
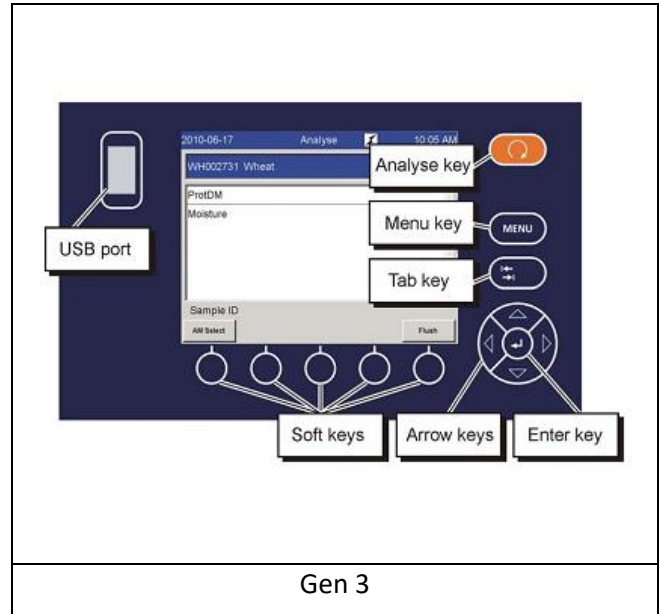


How to

Overview of Quick Menu

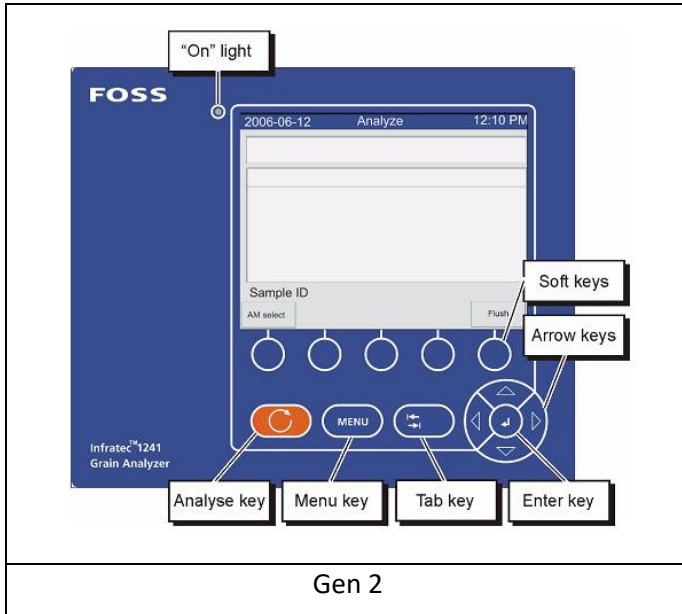

- 1 Main Menu
- 2 Instrument Information
- 3 Volume Information
- 4 Printer (On)
- 5 Test Weight Unit (On)
- 6 Save Scan (On)
- 7 Save Result (On)

Quick Menu Item	Function
Main Menu	Brings up User Menu (password protected)
Instrument Information	Shows the instruments serial number, software version, analyze counter, O- and P-constants and model.
Volume Information	Shows the amount of free space on the database partition.
Printer	Activates/deactivates the printer.
Test Weight Unit	Activates/deactivates the test weight module.
Save Scan	Stores scans for all AMs if activated. If deactivated, the AM scan storing settings will be used.
Save Result	Stores results for all AMs if activated. If deactivated, the AM result storing settings will be used.