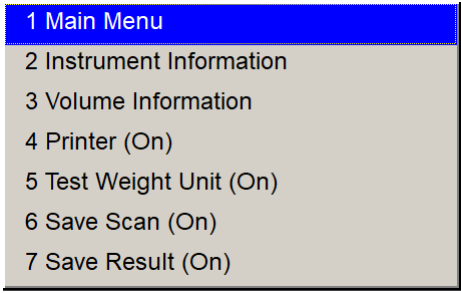
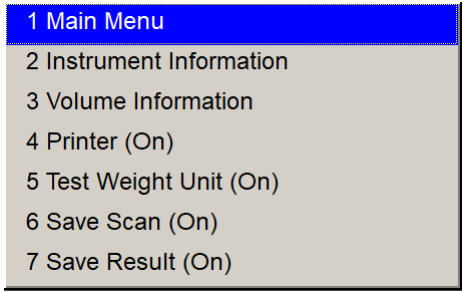
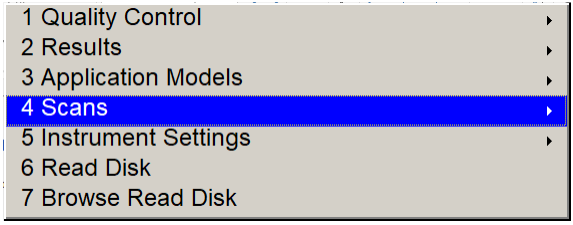
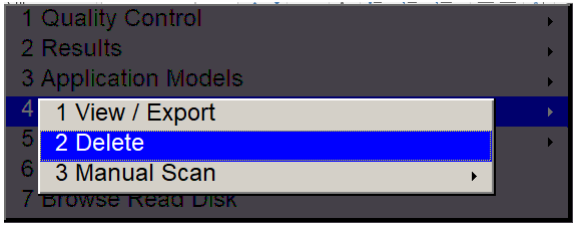
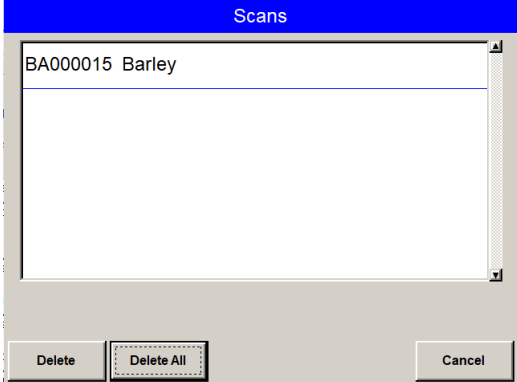
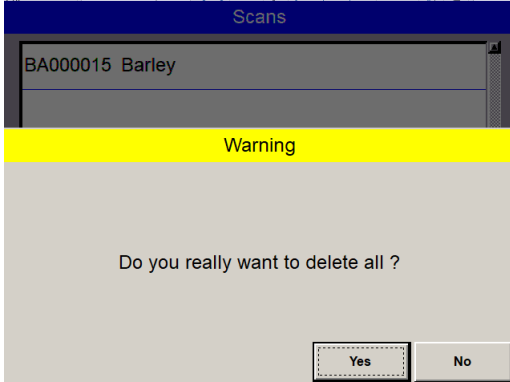
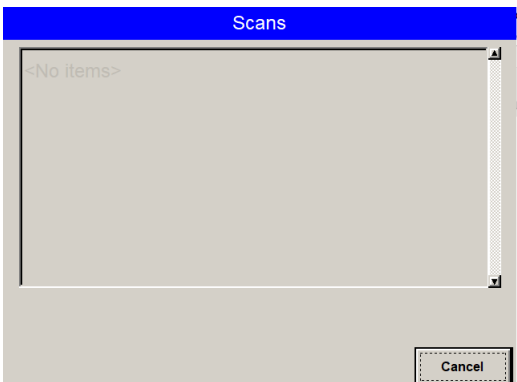
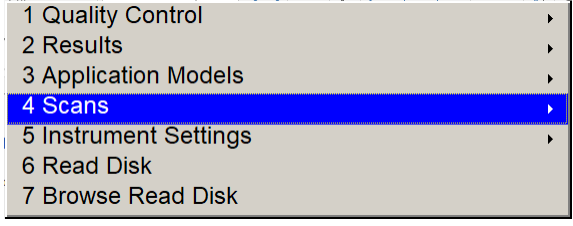









How to Delete Scans and/or Results

Before deleting scans/results, please synchronise to Mosaic manually to ensure all data is saved on the server

How to delete scans:

<p>1. Press MENU to access Quick Menu.</p> 	<p>2. Select "1 Main Menu". Enter password 123456.</p> 
<p>3. You are now in the Main Menu. Select "4 Scans"</p> 	<p>4. Select "2 Delete"</p> 
<p>5. Using the tab key, select "Delete All"</p> 	<p>6. Select "Yes"</p> 
<p>7. Press "Cancel"</p> 	<p>8. Press the left arrow to return to Main Menu. If you would like to delete results, see the next section. Otherwise, press MENU button to return to operator screen.</p> 

How to delete results:
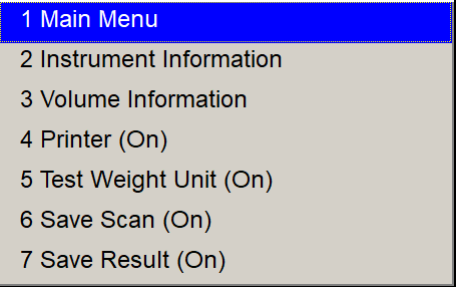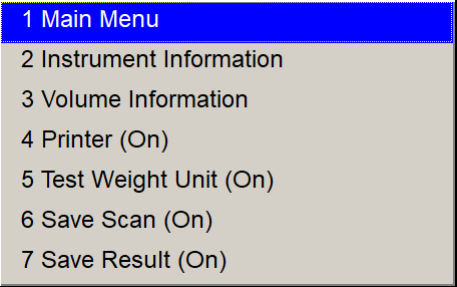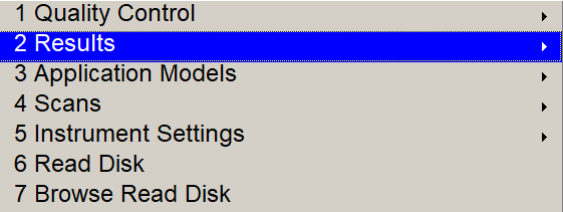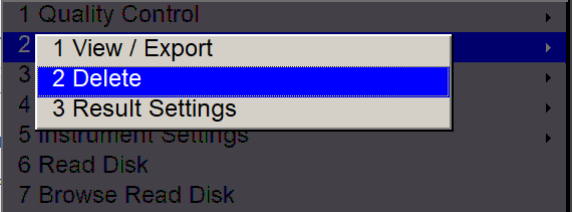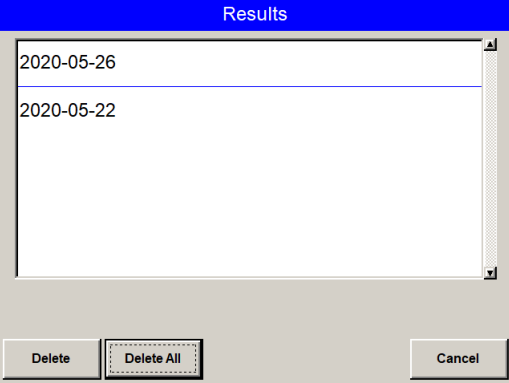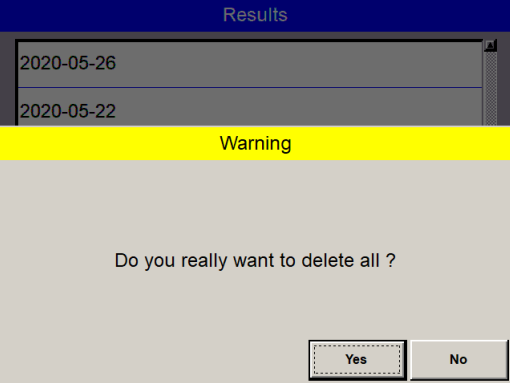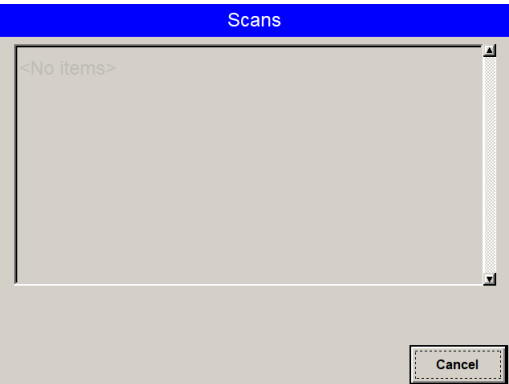
<p>1. Press MENU to access Quick Menu.</p> 	<p>2. Select "1 Main Menu". Enter password 123456.</p> 
<p>3. You are now in the Main Menu. Select "4 Scans"</p> 	<p>4. Select "2 Delete"</p> 
<p>5. Using the tab key, select "Delete All"</p> 	<p>6. Select "Yes"</p> 
<p>7. Press "Cancel"</p> 	<p>9. Press the left arrow to return to Main Menu. Press MENU button to return to operator screen.</p> 